


How to connect to LIS (Double communication)

2024.12.6 Made by After-sales Department


- 1、Click 'Setting--System--LIS', and setup LIS IP ' 192.168.2.110 ', Port ' 8080 ', Host IP ' 192.168.2.112 '

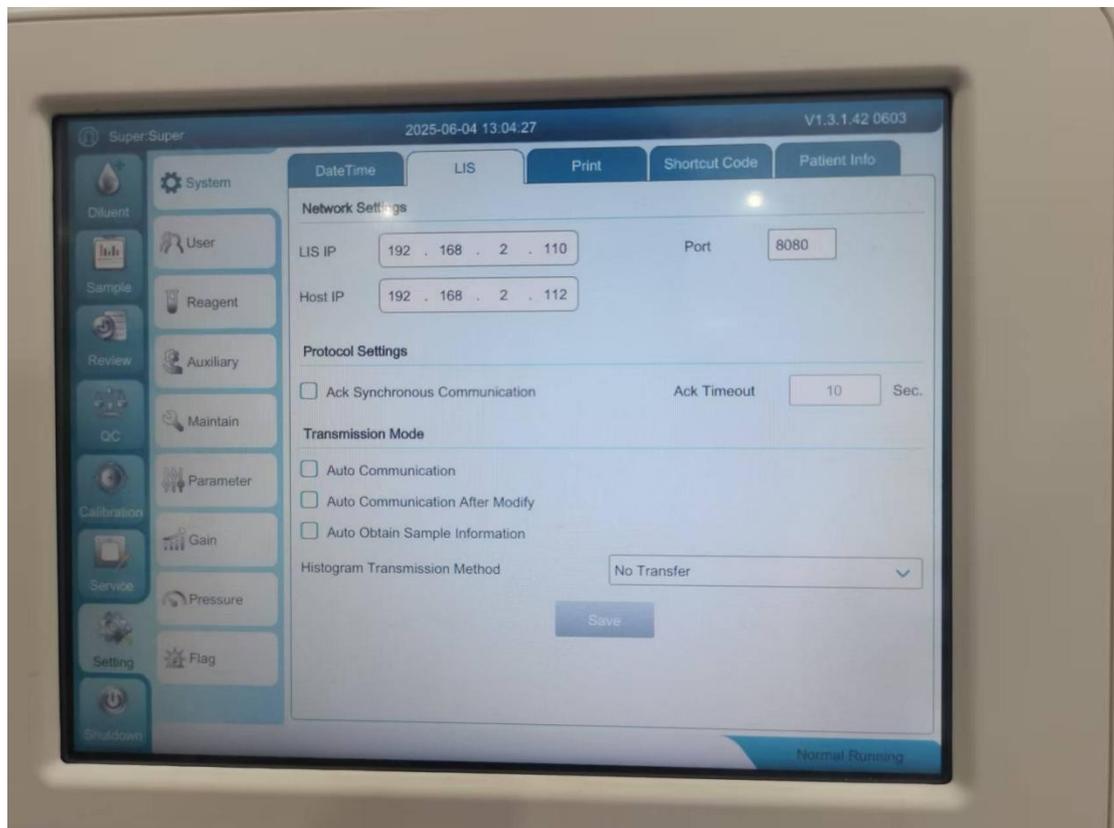


LIS IP is the IP of LIS system which you want to connect, Port is for communication

Host IP is the IP of the hematology analyzer

- 2、Open 'Hematology Data Management', click 'LIS Setup' and setup Hematology IP '192.168.2.112', Hematology '5100'

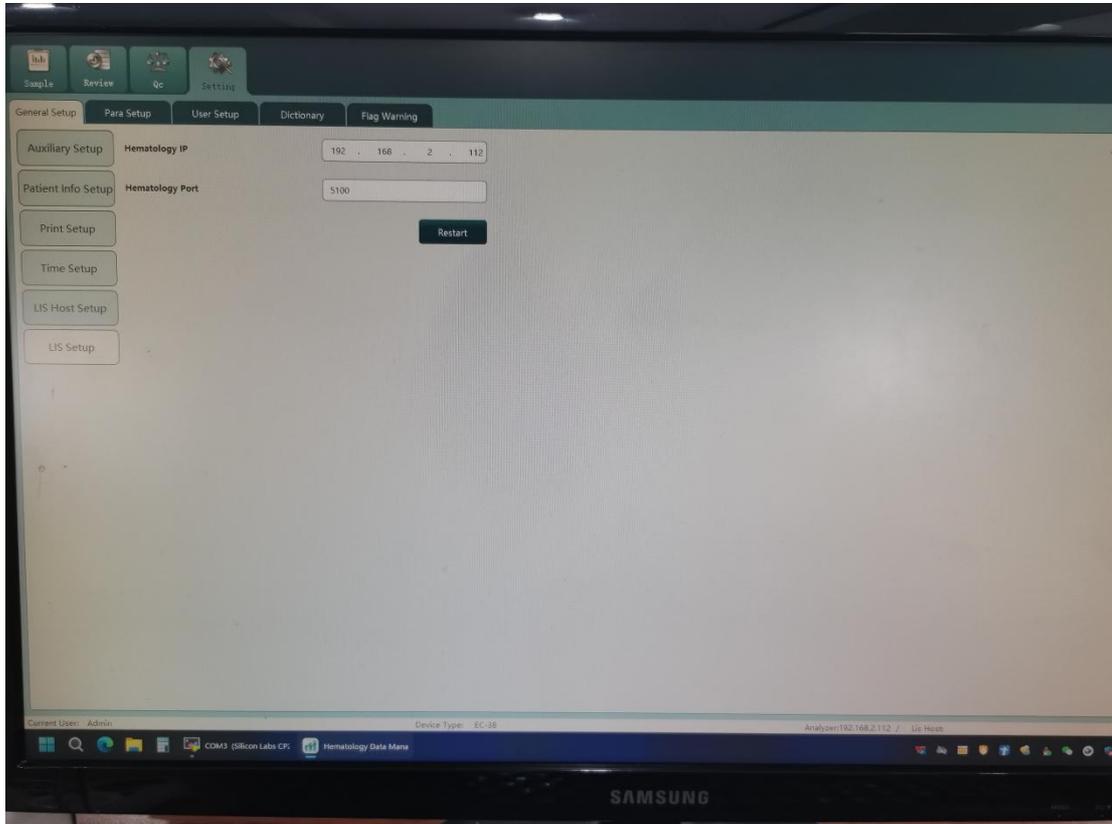

- 3、Click 'LIS Host Setup', check the Host Port whether is '8080'

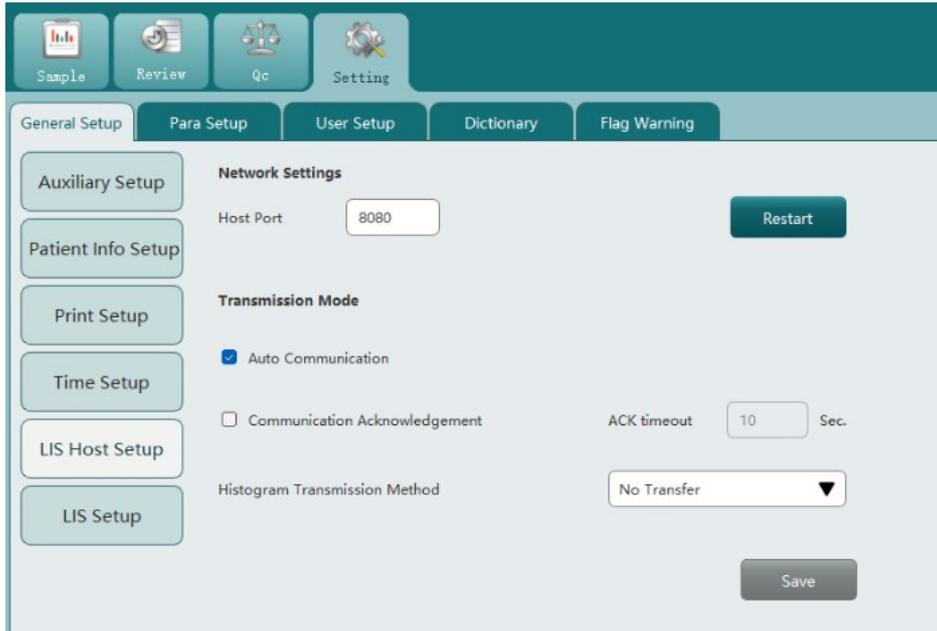

4、Set the computer IP setting, then connect the computer and the analyzer with a Ethernet cable.

